

Getting Started with the CBL2 and the DataMate Software

- Connect the CBL to a calculator
- Transfer DataMate to the calculator
 - Put the calculator in Receive mode, push TRANSFER
 - Start the DataMate Application
 - Press APPS, then ENTER, Main screen appears
- Connect the sensor(s)
 - Connect to correct channel
 - Calibrate the sensor
- Select the data collection mode.
 - On the main screen, press enter
 - Press the number beside the mode you want
 - Change the Time Graph Settings (optional)
 - Change the Advanced Time Graph Settings (optional)
- Collect the data
 - Press START
 - Store latest run - press TOOLS
- Graph the data
 - The Graph menu displays automatically
 - Press "up" or "down" arrow to move the cursor to the channel you want to view as a graph
- Analyze the data
 - On the DataMate main screen, press ANALYZE
 - Select the option you prefer
 - Press the number beside the option you want, then press ENTER
- Save, Retrieve, Delete Experiments
 - On the DataMate main screen, press ENTER, then Press SAVE/LOAD,
 - To Save: then SAVE EXPERIMENT, then key in the name, press ENTER
 - To Retrieve: then SAVE EXPERIMENT, then press the number you want, then ENTER
 - To Delete: then DELETE EXPERIMENT